# Innovate

Microbial Screening System

HARDWARE USER GUIDE

# contents

---

About Hygiena™ .....	1
Welcome to Hygiena's Innovate system .....	1
Document Conventions .....	1
Hygiena's Innovate system .....	1
<b>2.1</b> Safety Guidelines .....	1
<b>2.2</b> System Description and Technical Specifications .....	2
<b>2.2.1</b> Technical Specifications .....	2
<b>2.2.2</b> Plate Carrier .....	2
<b>2.2.3</b> Connections at Rear Panel .....	3
<b>2.2.4</b> Injectors .....	3
<b>2.2.5</b> Reagent Tray .....	3
<b>2.3</b> Innovate.im Requirements .....	3
Innovate system installation .....	4
<b>3.1</b> Space Requirements .....	4
<b>3.2</b> Connecting .....	4
<b>3.2.1</b> Remove the transport safety screw .....	4
<b>3.2.2</b> Electrical connections .....	5
<b>3.2.3</b> Reagent tray .....	5
<b>3.2.4</b> Waste and reagent tubing .....	5

## About Hygiena™

Hygiena™ is a world leader in the development and supply of rapid diagnostic and monitoring systems to detect and measure microbial contamination in finished products bound for consumers. The company has successfully developed tests for use in the dairy, beverage, personal care & cosmetic, and pharmaceutical markets and is associated with some of the best-known brands in the industry.

## Welcome to Hygiena's Innovate System

Welcome to Hygiena's Innovate system, designed to be the dairy and beverage industry standard for manufacturers using microbial screening to release their products to market. The Innovate system was developed based on input from dairy and beverage customers worldwide and is designed to offer simplicity of operation that is unmatched in the marketplace.

The Innovate system includes the Innovate bench-top device, Innovate.im software, RapiScreen™ reagent and service team. Together, all four components work together to provide the functionality and simplicity users need to meet their evolving demands.

## Document Conventions

The following conventions are used throughout this document to highlight specific information:

**Caution** – Warns the user that failure to avoid a specific action could result in physical harm to the user or the hardware.

**Caution** – Warns the user of a potential shock situation.

**Informational** – Advises of important information for completing the task.

## The Innovate System

Hygiena has done everything possible to guarantee that the equipment functions safely, both electrically and mechanically. The user must ensure the instrument is installed and maintained properly to guarantee safe operation.

The Innovate system has been manufactured in accordance with international safety requirements for electronic and medical measuring systems. The device is tested by the manufacturer and supplied in a condition that allows safe and reliable operation. Failure to follow the instructions may invalidate the warranty.

### 2.1 SAFETY GUIDELINES

Read this user's guide carefully before using the Innovate luminometer. Understanding the proper usage and operation is imperative for the safety of the user, as well as the protection of the instrument. The Innovate system luminometer must be installed and used in accordance with Hygiena recommendations. Installation must be performed by properly trained and authorized personnel.

Hygiena assumes no liability for any damages, including those to third parties, caused by improper use or handling of the instruments.

The instrument is manufactured in accordance to CE Marked, UL, CSA, EN 61010-1, EN 61326-1, EN 61000-3-2, and EN 61000-3-3 for electrical equipment.

It is the user's responsibility to have the instrument installed according to the local electrical codes (110 – 240V). The Innovate is equipped with a 3-wire grounded plug for safety. If your outlet does not accommodate the 3-wire plug, an electrician should install the correct outlet or use an adaptor to ground the instrument.

**DO NOT** open the instrument or tamper with any interior components. The instrument cover should never be opened while the instrument is running. Any replacement of parts shall be performed by the original manufacturer. Opening an instrument will void the warranty, and may cause a fire hazard or electric shock.

**DO NOT** place the instrument around any liquids, or use in a location where it may get wet. Liquids getting into the instrument could cause a fire hazard or electric shock. If liquid gets inside the instrument, unplug the power cord. Clean the unit or have it cleaned by an authorized service representative.

The system must only be used with Hygiena reagents and microtiter plates, in accordance with the instructions provided on the reagent kit insert.

Waste from the instrument's waste system (when priming/washing the instrument) must be disposed of properly. A plastic container is connected to the instrument to collect the waste. All statutory requirements for handling biological waste and reagents must be observed by the user.

The Innovate system instrument should be shipped in its own case. During transport the plate carrier must be secured by the transportation lock screw. The system needs to be washed and primed with washing solution leaving liquid in the injector lines. External tubing and injector needles should remain attached to the instrument and placed in a plastic bag.

The tests and maintenance work recommended by the manufacturer should be performed routinely to ensure the system operates as intended.

To obtain the best results, consider the following:

- Do not expose the instrument to direct sunlight.
- Install the instrument in a dry location where humidity levels do not exceed 85%.
- Keep plate carrier door closed at all times except when loading/unloading microtiter plates.
- Keep plate carrier free from dirt. Do not lean on plate carrier or move by hand.
- Clean up reagent spills on instrument immediately.
- Do not use the luminometer in an environment where the temperature exceeds 30 °C.

## 2.2 SYSTEM DESCRIPTION AND TECHNICAL SPECIFICATIONS

The Innovate system has been designed to deliver easy-to-use and robust capability to users. Due to its sophisticated technology and ease of use, it is well suited for dairy and beverage applications.

The instrument works with an extremely low-noise photomultiplier. The highest sensitivity possible is obtained due to the use of single photon counting technology and patented crosstalk reduction design. Specific Innovate system features include:

- Use of three (3) variable volume bellows-based injectors allowing precise, reliable, and automatic injection of reagent.
- Comprehensive shaking functionality to maximize mixing of sample and reagents.
- Temperature control allows reagents to be used up to five (5) days when stored in the Innovate system reagent cooling unit.
- Comprehensive Innovate.im software for instrument control and data management. Unique multi-tasking capability enables users to run one assay while preparing another.
- Connection to automated laboratory systems such as LIMS.
- Multiple user functionality allows multiple workstations to access results in the database via a network. Remote workstations can also access networked database results.

The Innovate system is a desktop microtiter plate luminometer and may be installed in any laboratory workplace meeting the manufacturer's recommended specifications.

### 2.2.1 Technical Specifications

Dimensions	33 x 43 x 27 cm
Operating Environment Conditions	Ambient temperature < 30 °C
Relative Humidity	< 85%
Voltage	110 – 240V
Weight	25.2 kg (Instrument, Reagent Tray, Cooling Unit)

### 2.2.2 Plate Carrier

The instrument contains a plate carrier accessible through the front panel door. Controlled via Innovate.im software, the plate carrier moves the microtiter plate to the defined measurement and injection positions.

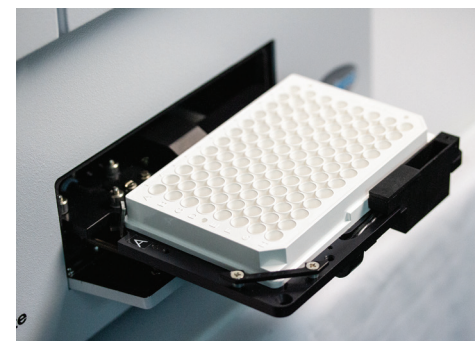
Figure 1: Plate carrier on instrumentation front panel.


### 2.2.3 Connections at Rear Panel

The PC connection, power supply, waste outlet, power switch, and instrument fuse are located on the instrument rear panel.

- USB port: Connects a computer to the Innovate system.
- Power plug: Connects the instrument to power using the cable supplied with the instrument.
- Power switch: Used to turn the instrument on and off.


Note: Turn the instrument on before starting the software, so the program can establish communication with the instrument.

- Waste outlet: Connects the waste container to the instrument.
- Fuse: Located next to the power switch in a black fuse holder.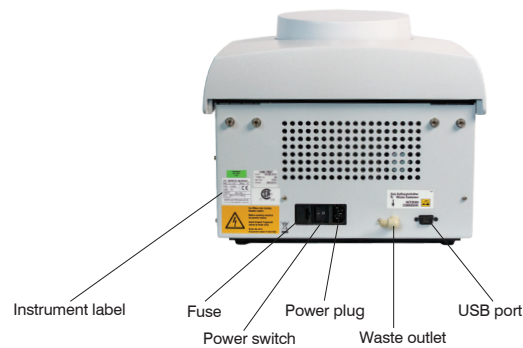

### 2.2.4 Injectors

The Innovate system has three variable injectors. Please see kit insert for setting up injectors with reagents. The injectors are located on the right side of the instrument. The tubing from the reagent vials is connected to the injector ports using screw-type caps.

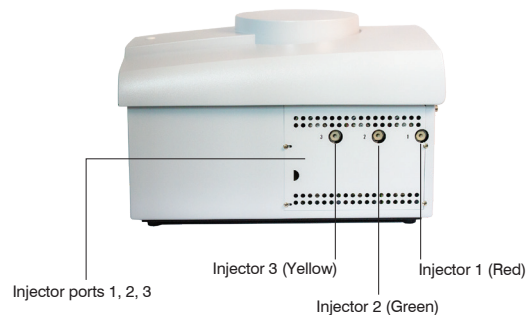### 2.2.5 Reagent Tray

The reagent tray is placed on the right side of the instrument. A special temperature control unit is placed inside the reagent tray in order to maintain optimal reagent temperature and prolong stability. Please see reagent kit insert for further details on reagent usage. Two washing solution vials are placed in the reagent tray to wash the injectors daily during startup and shutdown.

## 2.3 Innovate.IM SOFTWARE

### 2.3.1 Innovate.im Requirements

Innovate.im requires the following minimum computer hardware specifications:

Processor	1 GHz
RAM	1 GB
Hard Drive	100 MB free space
Graphics Card Resolution Support	1024x768
Optical Drive	CD-ROM
Operating System	Windows 7 Professional, Windows XP Professional, Windows Vista Business
Communication Ports	USB, Parallel (if needed for printer)
Web Browser	Internet Explorer 6.0 or higher (optional)
Peripherals	Mouse, Keyboard, Monitor

Multuser functionality requires the following minimum specifications in addition to the above system requirements. All users must have full read and write permissions for the networked folder where the database tables are stored. For optimal speed, no more than four systems should access the networked database tables at one time.

Network Speed	1 Gbps
RAM	2 GB
License	Needed for remote users not attached to an Innovate luminometer

IMPORTANT: Plate preparation files, report templates, pre-defined queries, and database export drivers must be configured identically on all networked systems.

## Innovate System Installation

This section will lead you step-by-step through the Innovate system luminometer installation process. The Innovate system should be set up in a dry, dust-free room and protected from direct sunlight and significant temperature fluctuations. It should not be set up next to heating or air conditioning units. The Innovate system luminometer should be operated in an environment where the temperature does not exceed 30 °C and humidity levels are below 85%.

### 3.1 SPACE REQUIREMENTS

The luminometer has the following dimensions:

- W x D x H: 33 x 43 x 27 cm
- The Innovate system luminometer should be set up so the rear panel connection ports are easily accessible and not subjected to physical stress. Allow for sufficient space on the side of the instrument for the reagent tray and the PC used for Innovate.im.

The original packaging is reusable and should be retained for any future transport of the instrument. When unpacking the instrument, make sure the shipment is complete and shows no sign of transport damage. The packaging is designed to eliminate transport damage. Should the instrument or instrument parts appear damaged, please notify your local Hygiena representative immediately.

Please verify that all components were shipped and received in good condition by referencing the packing list provided with the instrument. If items are missing or damaged, please contact Hygiena immediately.



## 3.2 CONNECTING

### 3.2.1 Remove the transport safety screw

The red label is placed in front of the Innovate system luminometer to remind you to remove the transport safety screw (a). Remove the white cap and unscrew the instrument cover with the supplied screwdriver (b); unlock the lid by loosening the screw. Open the lid.

Remove all of the foam pieces inside the instrument (c, d).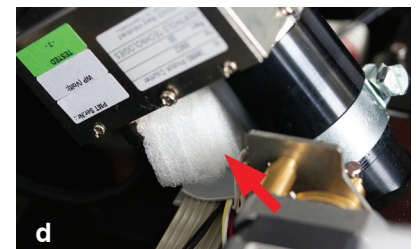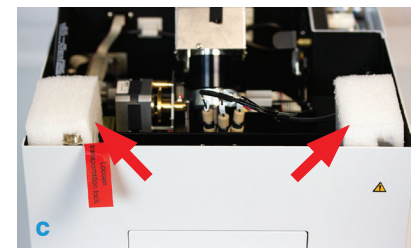

Unscrew the transport safety screw (e). This screw secures the plate carrier and should always be engaged to prevent damage when shipping the instrument.

Place the red label inside the instrument (f) to the right side of the transport screw.

Close the instrument cover and replace lid screw and white cap (g).


### 3.2.2 Electrical connections

Connect the USB cable provided to the port located on the rear of the Innovate system. Do not connect the USB cable to the PC until installation of Innovate.im software. (See Section 3.1).

Note: Check if power supply is within the permissible range of the Innovate operating voltage. Connect instrument only if the power supply is 110 – 240 V.


- Connect instrument to the wall outlet using the power cord supplied.
- If you have not already done so, connect the computer to the wall outlet.
- Connect temperature control unit to the wall outlet using the power cord supplied.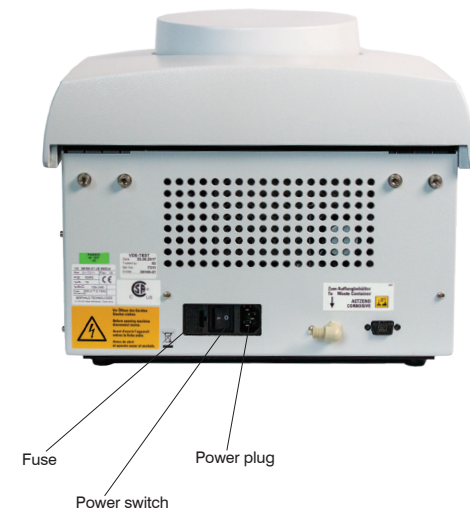

### 3.2.3 Reagent tray

Assemble the reagent tray using the Allen key supplied with the Innovate system. The tray contains two tube holders, one with an opening and one without an opening. Place the tube holder with the opening at the front side of the tray.

The liquid level will be visible through the opening. Place the cooling unit inside the tray.

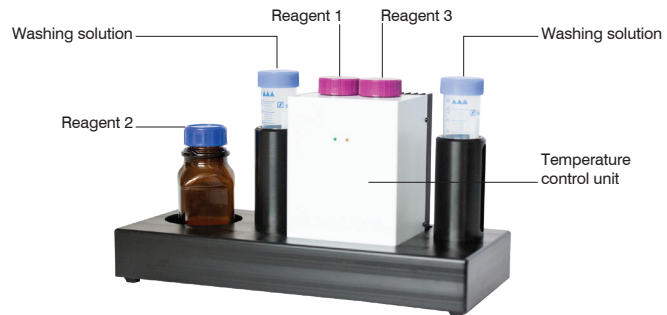
### 3.2.4 Waste and reagent tubing

Connect supplied waste tubing to the waste outlet.

- Add a few drops of antifoam to the waste container.
- Place waste tubing into the waste container to collect the waste solution.
- Instrument is shipped with external reagent tubing connected to injectors 1, 2, and 3.

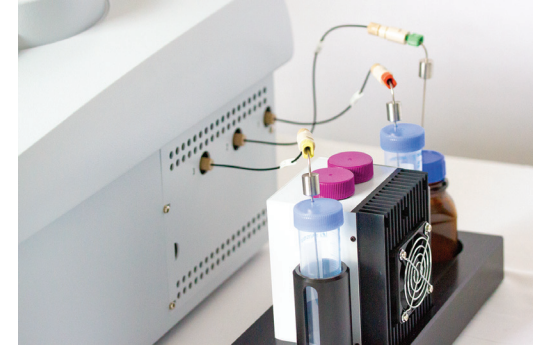- Place all three injector needles in washing solution.
- Note: Wear gloves when touching the injector needles.

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